




***Midwest LRIG Rapid Fire***  
Solutions for Automated Liquid Handler  
Optimization

*April 21, 2022*

*Dana Campbell*  
*Senior Applications Specialist*

# Global Standard for Liquid Handling Performance

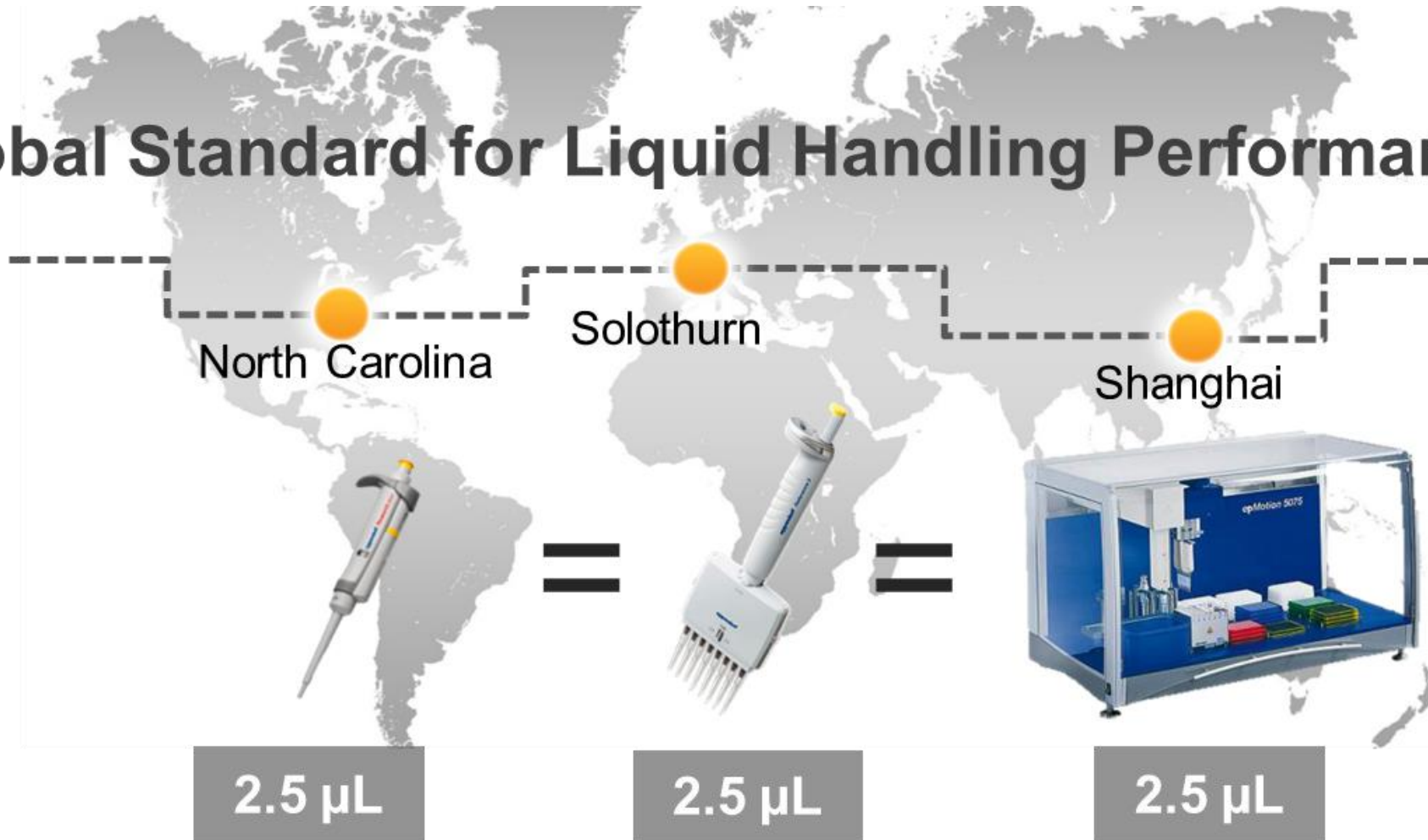

# MVS – Multichannel Verification System

*Calibration, Verification, Operator Training and Liquid Class Optimization*


- To help you with your ALH optimization, there are 4 different QualAssure solutions, so you can choose the ones that best mimic your liquid types:

- Aqueous

- DMSO


- PCRMix


- SerumSub

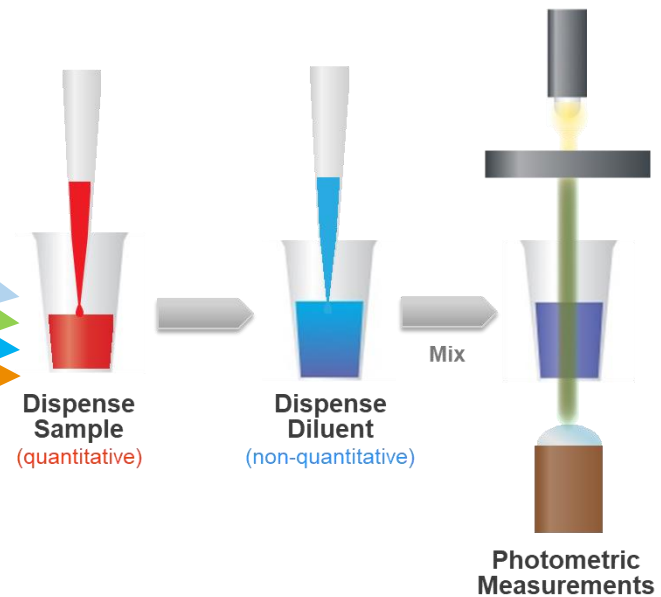

# Why Perform Liquid Class Optimization

*Reagents and Samples are Characterized by their physical properties (liquid class parameters)*

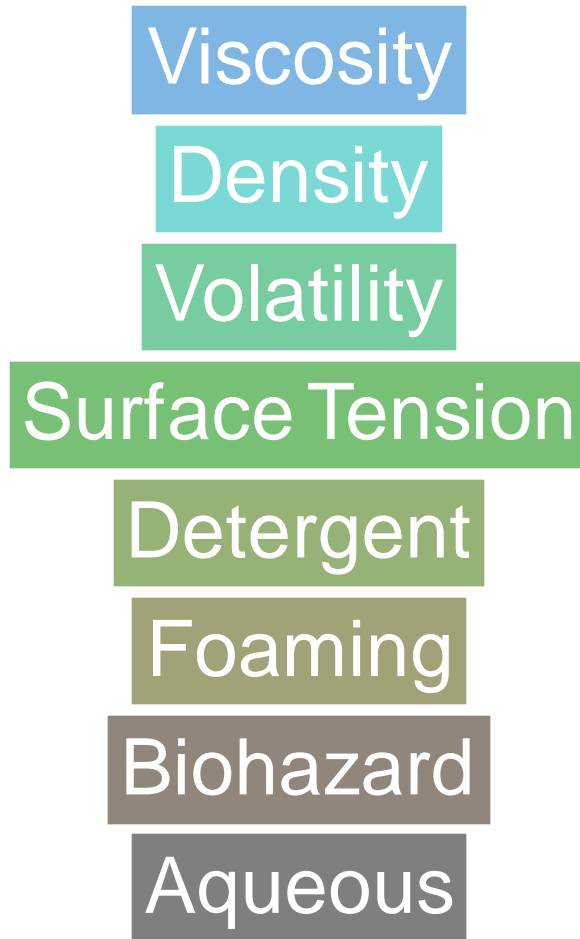

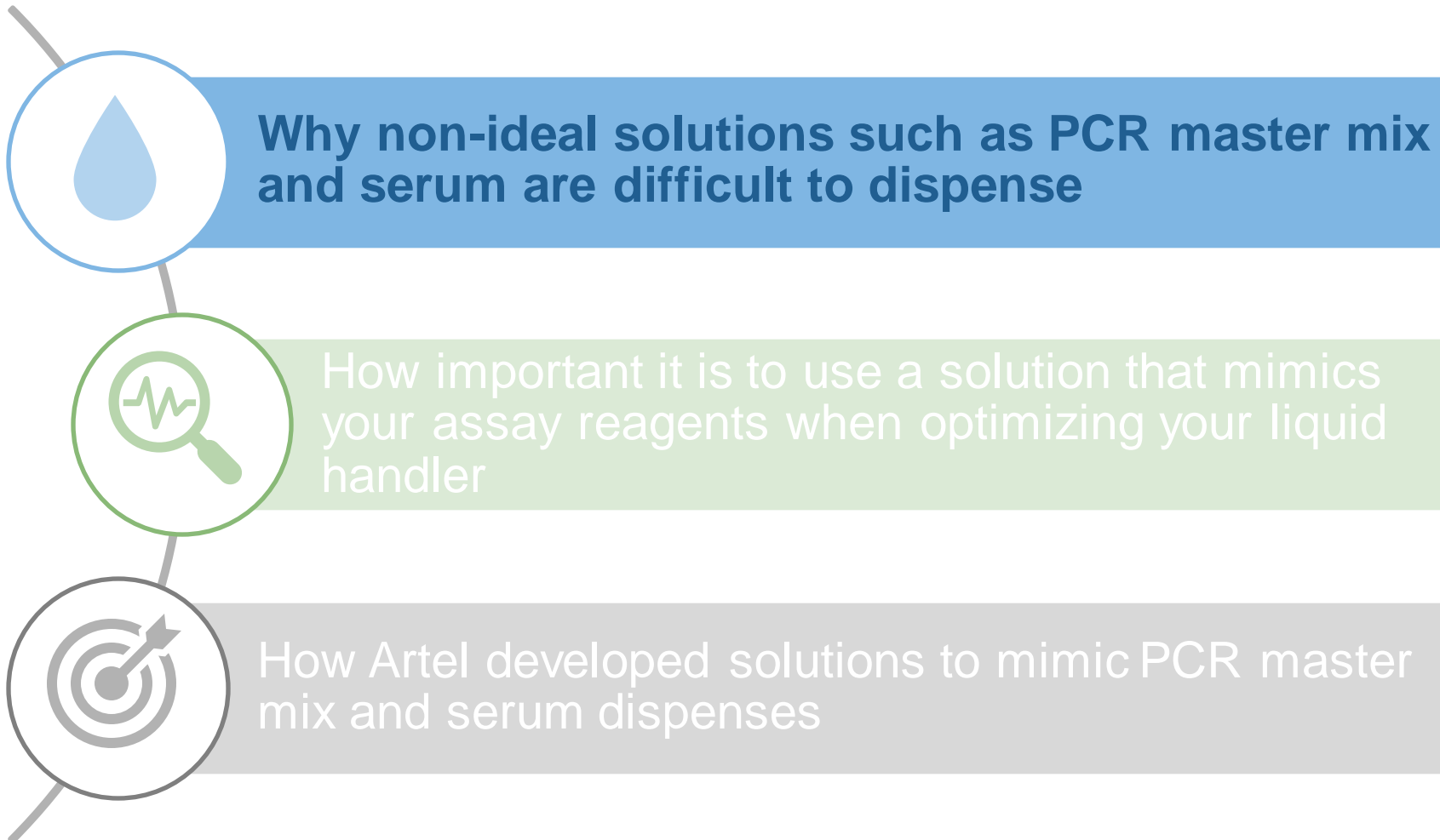

**Why non-ideal solutions such as PCR master mix and serum are difficult to dispense**

How important it is to use a solution that mimics your assay reagents when optimizing your liquid handler

How Artel developed solutions to mimic PCR master mix and serum dispenses

# Not all liquids behave the same

*Many parameters affect pipettability*


- Viscous
- Dense
- Vapor Pressure
- Volatile
- Detergent
- Volume dependent
- Foaming
- Surface Tension



# What is serum?

- What is it?


- The fluid obtained after the complete coagulation of the blood, without adding an anticoagulating agent.

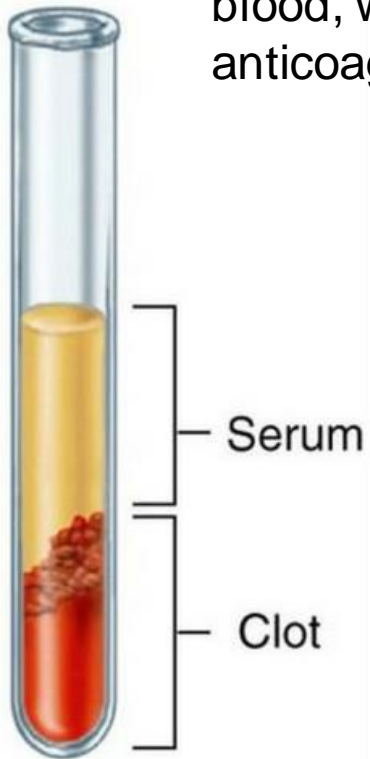

- What's in it?

- Glucose → Increases viscosity
- Hemoglobin
- Albumin (~50%) → Protein, reduces surface tension
- Immunoglobulins (~35%)
- Cholesterol → Increases stickiness
- Triglycerides
- Sodium
- Calcium
- Chloride
- Potassium
- And other components

High protein content increases bubbles and foaming

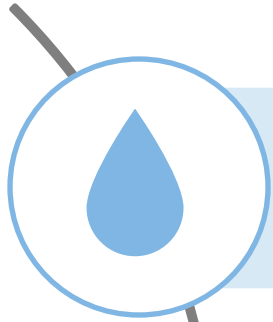

Why non-ideal solutions such as PCR master mix and serum are difficult to dispense




**How important it is to use a solution that mimics your assay reagents when optimizing your liquid handler**



How Artel developed solutions to mimic PCR master mix and serum dispenses



# Poor Accuracy and Precision Impacts Assay Results

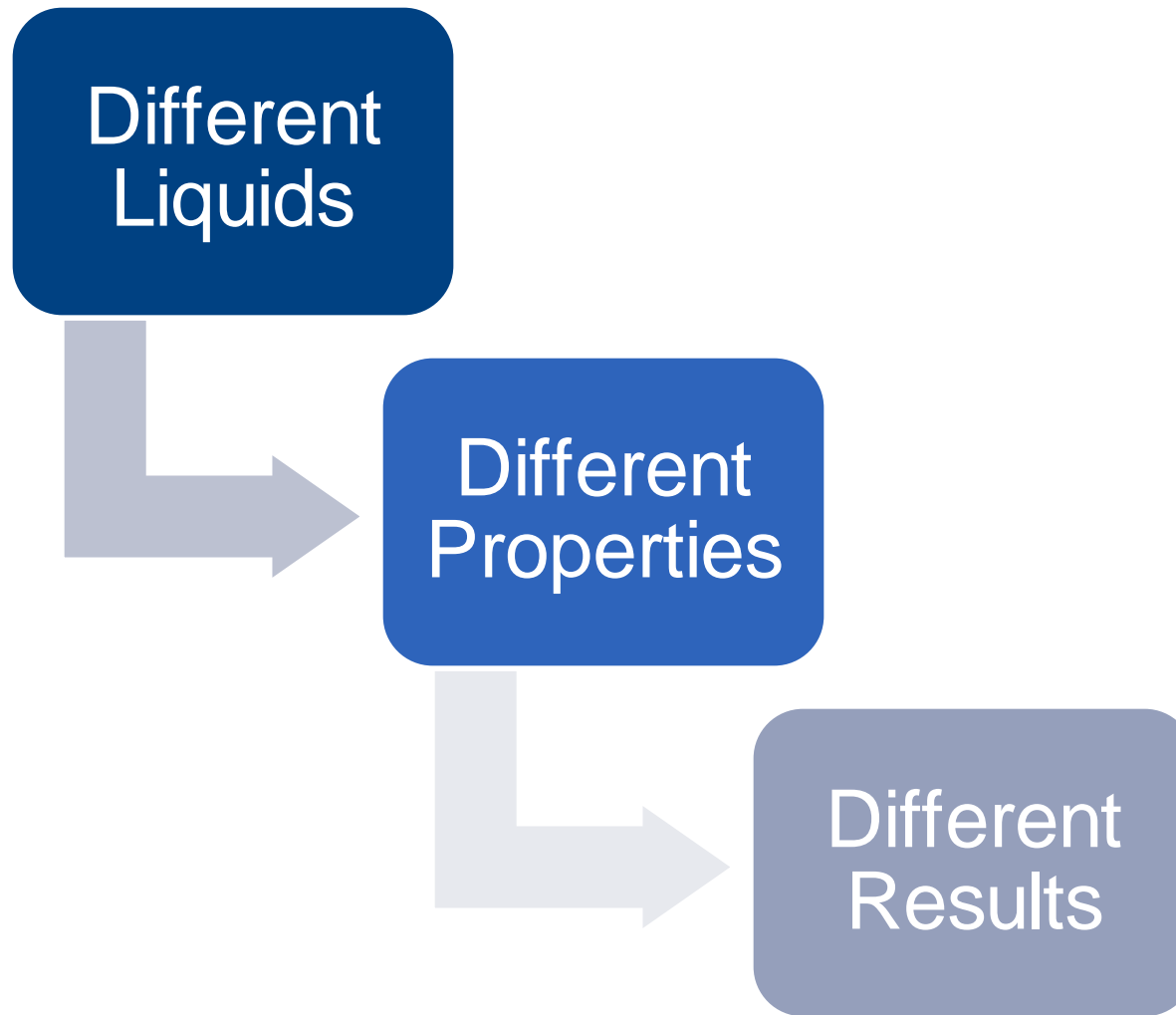

# Benefits of Using SerumSub

---

- Most clinical samples require an extraction step and dilution step, so quantitative transfer is essential
- Limited availability of serum sample
- Increase the accuracy and precision of liquid transfers
- Speed up and improve development and transfer of assays
- Minimize false negatives and false positives
- Minimize overall assay variability
- Maximize data quality and reproducibility
- Ability to test assays daily.
- Improve efficiency by reducing assay transfer problems
- Liquid class optimization for automated liquid handlers

# Process Optimization

*Not all liquids are created equal*



Liquid Class Optimization



Liquid Handler Evaluation



Calibration



Operational & Performance Qualification



Operator Training



Method Transfer & Scale Up



Interim Verification

## What is Liquid class optimization?



- Adjustment of any parameters on an automated liquid handler are intended to:
  - **increase** the accuracy and precision of liquid transfers
  - **decrease** variability of assay results and even false results.


# Liquid Class Optimization

Before and After...




# Aqueous Liquids vs Serum “like” Liquids

*Serum and Aqueous liquids do not behave the same*

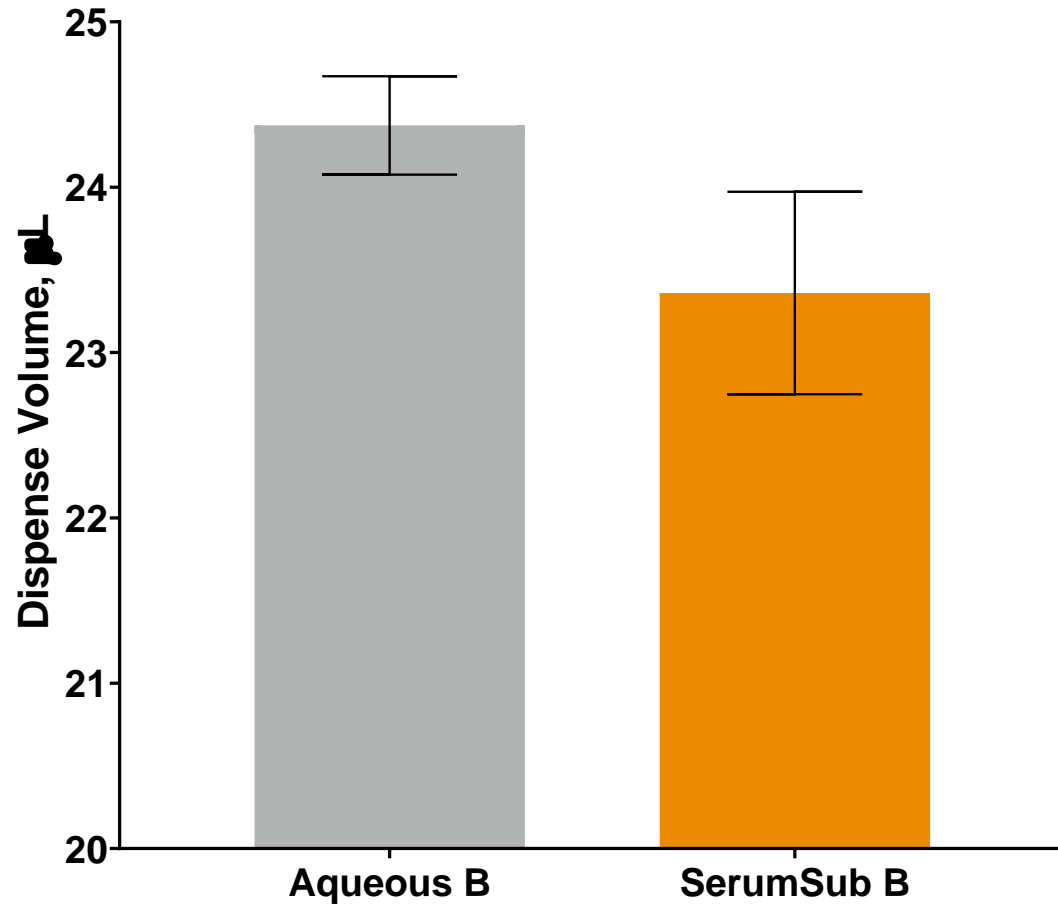


# Liquid Class Optimization – Simplified

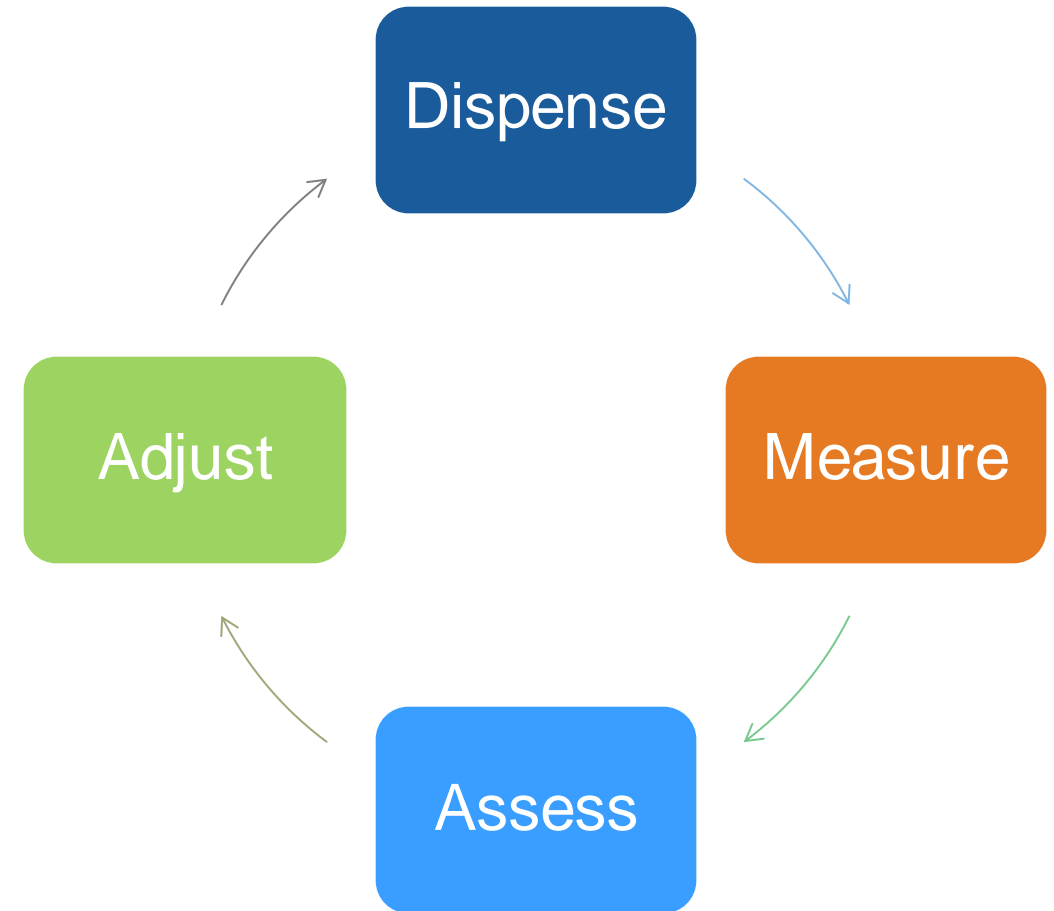


# Liquid Class Recommendations for SerumSub

Start Here

- Use a slower aspiration & dispense speed
- Add or increase a pause/settling time
- Add tip touch
- Continue adjusting 1-2 parameters at a time until ideal settings are determined

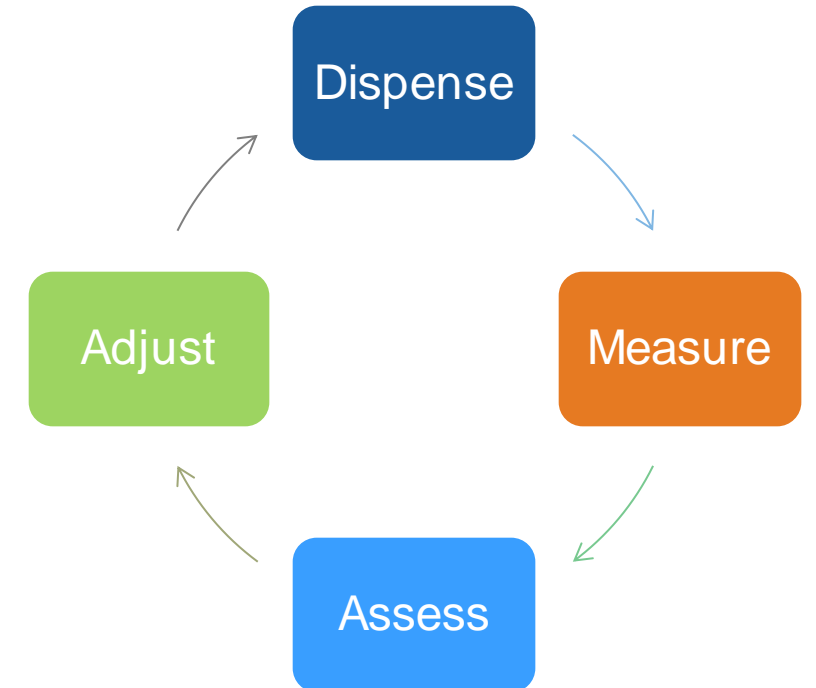



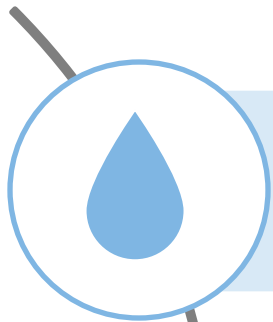

Why non-ideal solutions such as PCR master mix and serum are difficult to dispense



How important it is to use a solution that mimics your assay reagents when optimizing your liquid handler




**How Artel developed solutions to mimic PCR master mix and serum dispenses**

# Equivalency Testing

A test to determine if one is like the other.

# Equivalency Testing

A test to determine if one is like the other.

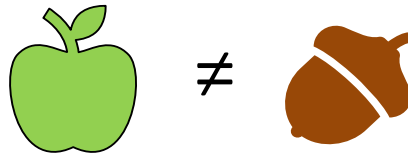




# Equivalency Testing

A test to determine if one is like the other.






# Equivalency Testing

A test to determine if one is like the other.

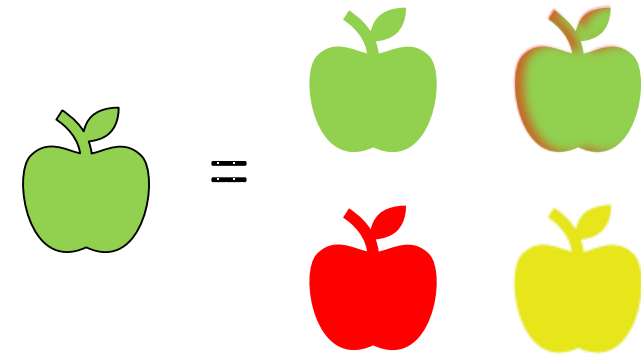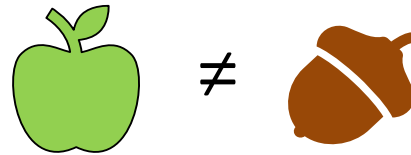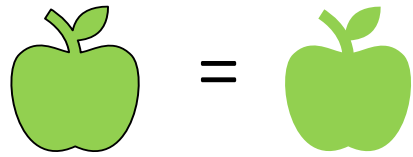

# Equivalency Testing




A test to determine if one is like the other.

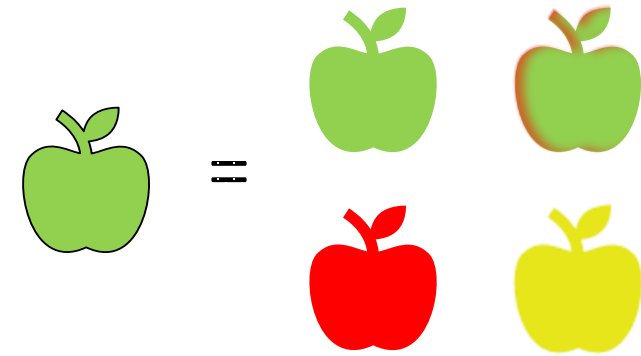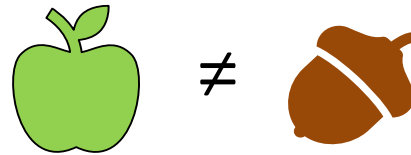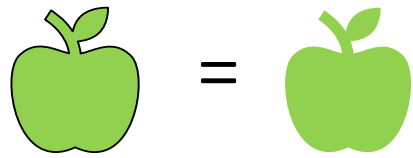

1. Identify different master mixes or sera to test.

# Equivalency Testing




A test to determine if one is like the other.

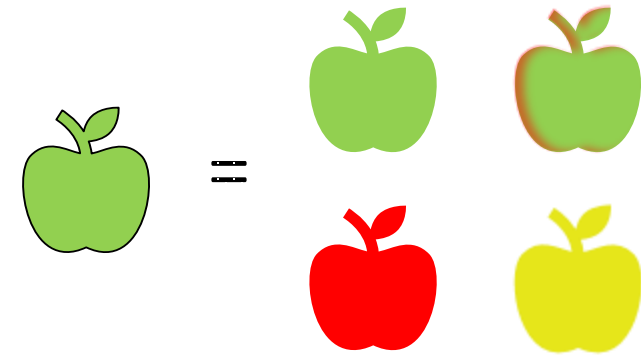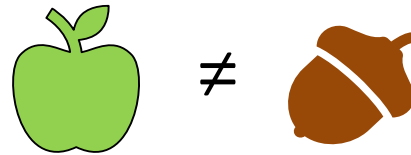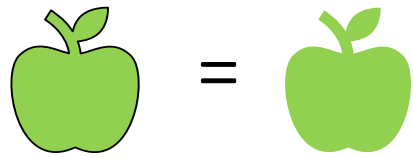

1. Identify different master mixes or sera to test.
2. Develop simple transfer method on an ALH.

# Equivalency Testing


A test to determine if one is like the other.

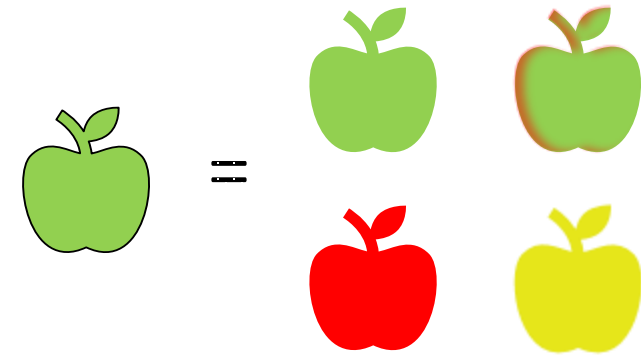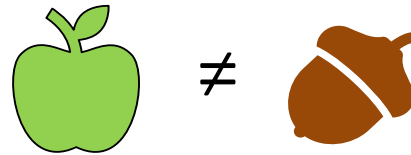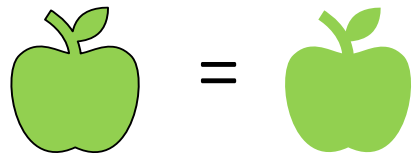


1. Identify different master mixes or sera to test.
2. Develop simple transfer method on an ALH.
3. Use gravimetry to measure and compare.



## Comparison to Commercial Sera

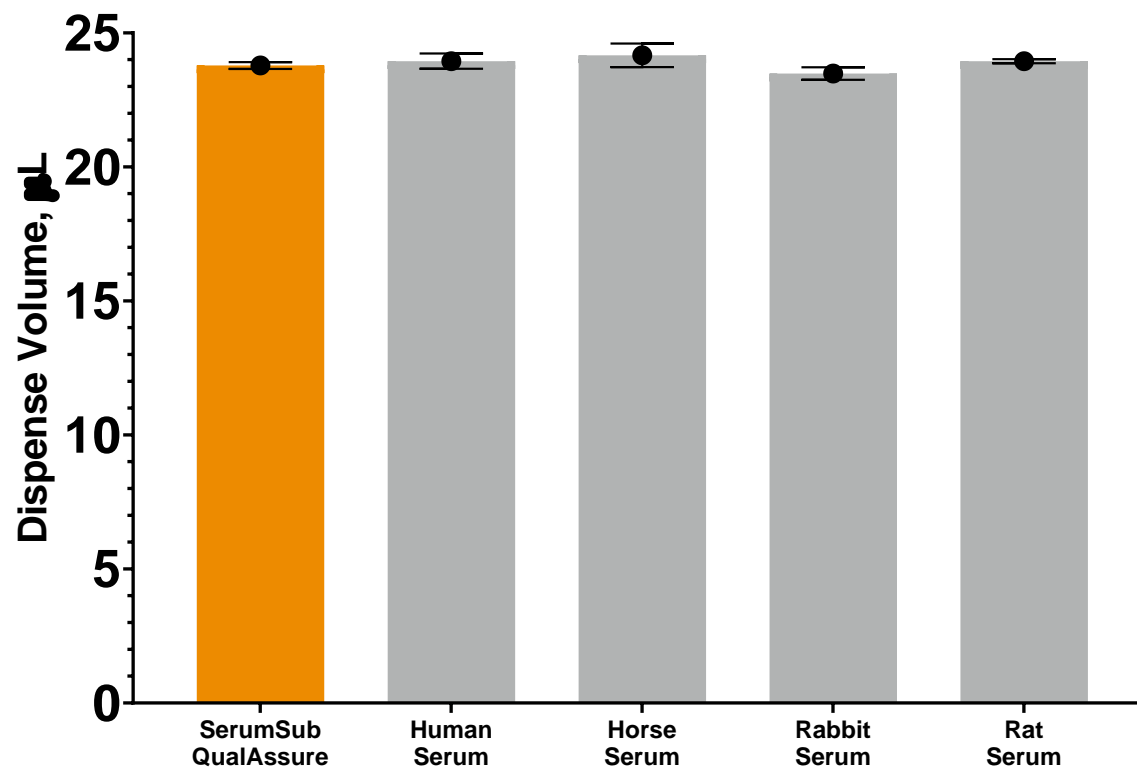

- SerumSub mimics dispensing performance of various animal sera
- Allows users to simulate and optimize dispensing steps without wasting expensive reagents

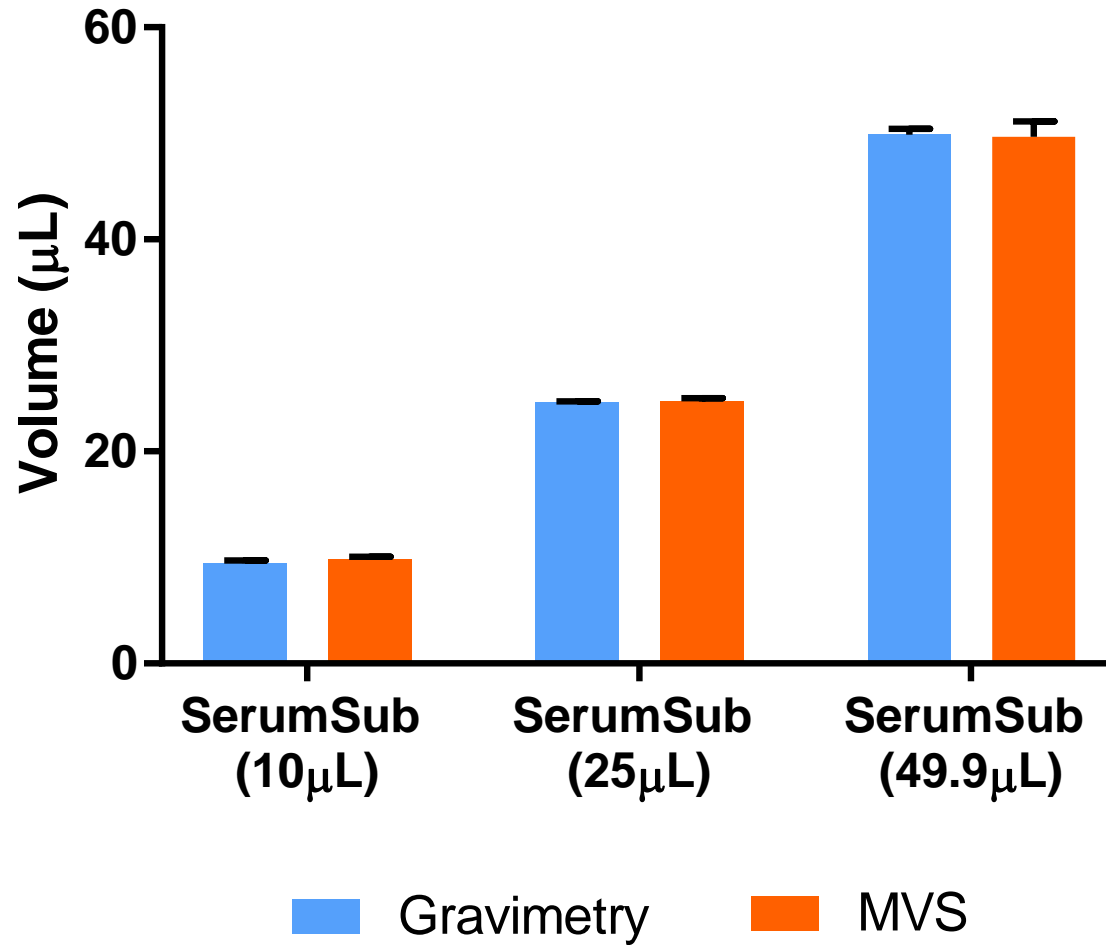



PCRMix


2 to 49.99  $\mu$ L in 96-well plates

SerumSub

10 to 200  $\mu$ L in 96-well plates



# Summary

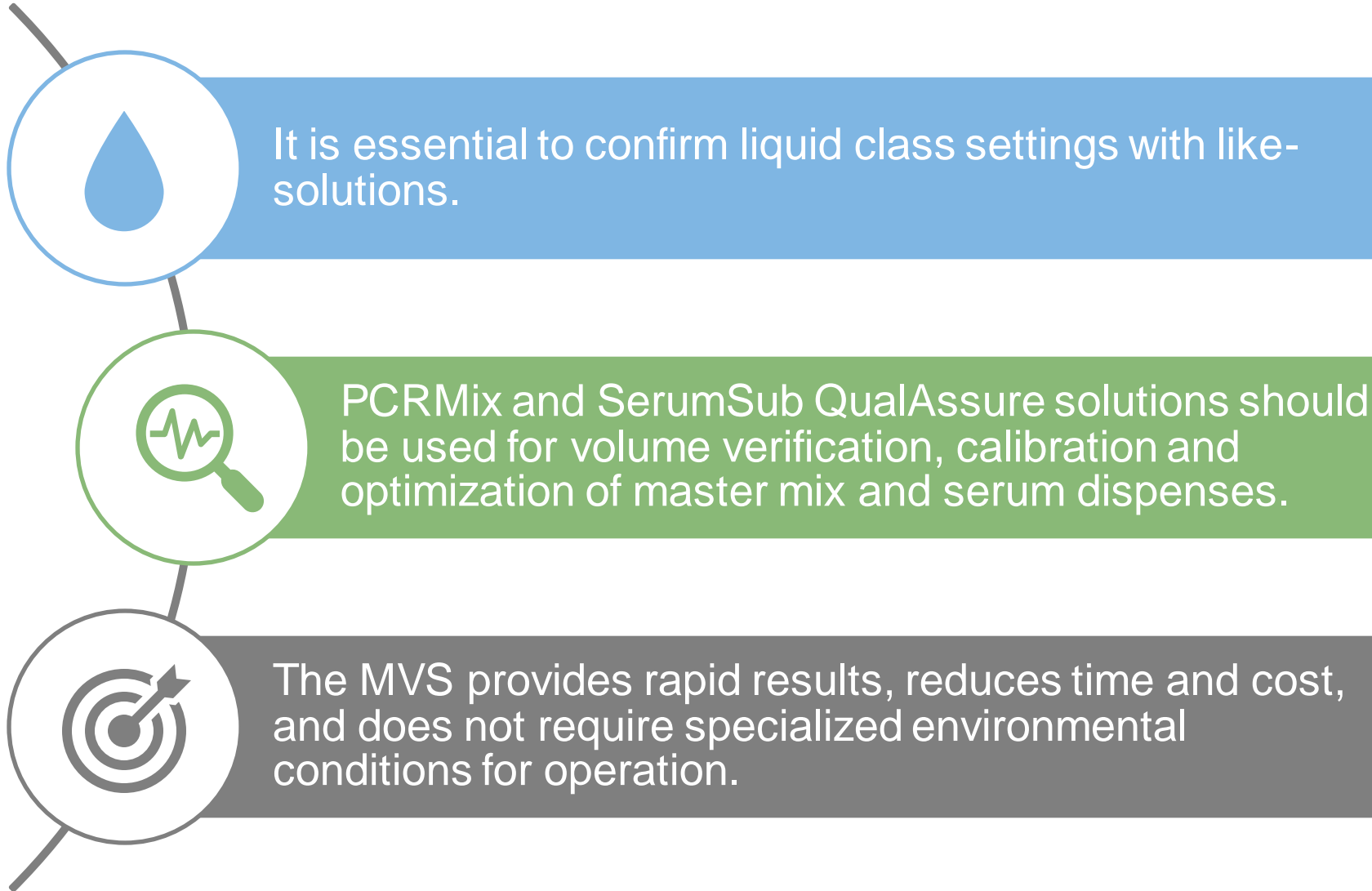



Thank you very much for your attention!

Contact me at: [dcampbell@artel.co](mailto:dcampbell@artel.co)