

## MVS Data Manager Release Notes v3.7.0.2

### Feature Changes

- Added functionality for PlasProxy QualAssure™ A and B in 96-well plates
- Added MVS System Specifications for PlasProxy:

#### MVS System Specifications for PlasProxy A and B in 96-well Verification Plates

Volume Range	Inaccuracy	Imprecision
10.00 µL to 200.0 µL	5.0%	0.80%

### Compatibility Changes

- Data Manager 3.7.0.2 is required for use of PlasProxy.
- PlasProxy is not supported in 384-well Standard Profile, 384-well Low Profile, or 384-well Low Volume plates.
- An update to Large Volume Calculator version 2.10 is required to create summations with DM 3.7 verification reports.

### Bug Fixes

There are no required bug fixes with this version.

### Updating

Please refer to the **MVS System Upgrade 3.7 – IQOQ** document for instructions to update Data Manager software.

If you have any additional questions or concerns, please contact Artel Technical Support at 888-406-3463 option 2 or by email at [support@artel.co](mailto:support@artel.co).

7A7368A